

US ENVIRONMENTAL PROTECTION AGENCY OFFICE OF PESTICIDES PROGRAMS REGISTRATION DIVISION (TS-767) WASHINGTON, DC 20460	EPA REGISTRATION NO. 46781-5	DATE OF ISSUANCE NOV 18 1988
	TERM OF ISSUANCE	
	NAME OF PESTICIDE PRODUCT ... Machine Concentrate	

NOTICE OF PESTICIDE:  REGISTRATION  
 REREGISTRATION  
 (Under the Federal Insecticide, Fungicide and Rodenticide Act, as amended)

NAME AND ADDRESS OF REGISTRANT (Include ZIP code)

NOTE: Changes in labeling formula differing in substance from that accepted in connection with this registration must be submitted to and accepted by the Registration Division prior to use of the label in commerce. In any correspondence on this product always refer to the above U.S. EPA registration number.

On the basis of information furnished by the registrant, the above named pesticide is hereby Registered/Reregistered under the Federal Insecticide, Fungicide, and Rodenticide Act.

A copy of the labeling accepted in connection with this Registration/Reregistration is returned herewith.

Registration is in no way to be construed as an indorsement or approval of this product by this Agency. In order to protect health and the environment, the Administrator, on his motion, may at any time suspend or cancel the registration of a pesticide in accordance with the Act. The acceptance of any name in connection with the registration of a product under this Act is not to be construed as giving the registrant a right to exclusive use of the name or to its use if it has been covered by others.

The following conditions apply to the registration of this product with the Agency (Section 6(a) of the Act):

1. Since you are a manufacturer of a pesticide registered under Federal Insecticide, Fungicide, and Rodenticide Act, you are required to register any similar products to be sold in interstate commerce.

2. Use the phrase "EPA Registration No. 46781-5" on your label when you release the product for shipment.

3. Submit three (3) copies of your final printed labeling before you release the product for shipment, with the enclosed enclosure for a further description of final printed labeling.

If these conditions are not complied with, the registration will be subject to cancellation in accordance with 40 CFR section 6(e). Your release for shipment of the product constitutes acceptance of these conditions.

A stamped copy of the label is enclosed for your records.

John H. Lee  
 Product Manager (31)  
 Antimicrobials Programs Branch  
 Registration Division (TS-767C)

**BEST AVAILABLE COPY**

enclosure

ATTACHMENT IS APPLICABLE

SIGNATURE OF APPROVING OFFICIAL

DATE

METR

# AUTOMATIC MAC

## A disinfectant for use in auto

For complete disinfection of medical equipment, thoroughly clean, rinse, and rough-dry objects before immersing in METRICIDE Automatic Machine Concentrate. Cleanse and rinse the lumens of hollow instruments before using with METRICIDE Automatic Machine Concentrate.

**Disinfection:** Immerse instruments or equipment completely for a minimum of 10 minutes at 20°C. to destroy vegetative organisms including *Pseudomonas aeruginosa*, *Staphylococcus aureus*, *Salmonella choleraesuis*, and pathogenic fungi as represented by *Trichophyton mentagrophytes*. For 100% kill of *Mycobacterium tuberculosis*, immerse instruments or equipment completely for a minimum of 30 minutes at 30°C., 60 minutes at 25°C. or 70 minutes at 20°C. **VIRUSES:**

The following viruses are completely inactivated in 3 minutes at 20°C.: *Respiratory Syncytial Virus*; *Adenovirus 2*; *Coxsackievirus B5a*; *Echovirus 8*; *Poliovirus 1 & 2*; *Vaccinia virus*; *Influenza A<sub>1</sub> HK*. After immersion, remove instruments or equipment from METRICIDE Automatic Machine Concentrate and rinse thoroughly. All times and temperatures are based on EPA-approved testing on inanimate surfaces.

1. Dilute contents of three (3) 1-gallon containers of METRICIDE Automatic Machine Concentrate with twelve (12) gallons of cold tap water.
2. Add contents of half-gallon container of Activator to solution.
3. **Act.** When activated, METRICIDE Automatic Machine Concentrate changes color to yellow and is ready to use.

Product No. 4  
**MX-1480**  
 Active Ingredient  
 Glutaraldehyde ..... 12.5%  
 Inert Ingredients ..... 87.5%  
 Total ..... 100.0%  
 Contains Sodium Nitrite as a corrosion inhibitor.

1. Follow directions for use of your Automatic Decontamination Machine.
  2. Do not use until directions for activation have been followed.
- Do not use activated solution beyond 14 days.  
 It is a violation of Federal Law to use this product in a manner inconsistent with its labeling.

**DANGER: Keep out of reach**  
 See right side panel for precautions  
 Note: Activating solution must be added  
 See directions for activation

E.P.A. Reg. No. 46781-2 L  
 E.P.A. Est. No. 46781-CO-1

ACCEPTED  
 with COMMENTS  
 in EPA Letter Dated:

NOV 10 1988

Under the Federal Insecticide,  
 Fungicide, and Rodenticide Act as  
 amended, for the pesticide  
 registered under EPA Reg. No.

46781-5

**BEST AVAILABLE COPY**

3/3  
**ICIDE™**

# LINE CONCENTRATE

ting solution  
omatic machines

of children.  
utions.

to concentrate before this product is effective.  
\*See directions for use

3  
**DANGER: Hazard to Humans and Domestic Animals**  
Follow All Label Directions

Eyes - Safety glasses or goggles are recommended. Contact with eyes causes damage. In case of contact, flush with water immediately and get medical attention.

Skin - Avoid skin contact as possibility of sensitization exists. May cause skin irritation. Use gloves. For skin contact flush immediately and thoroughly with water.

Ingestion - Avoid contamination of food. Harmful if swallowed.

Inhalation - Use in ventilated area.

Keep in a cool place. Do not reuse empty container. Wrap container and put in trash.

Especially recommended for disinfection of . . .

Anesthesia equipment  
Respiratory therapy equipment  
Metallic equipment or instruments

Rubber objects  
Plastic objects  
Thermometers

Contents One Gal. (128 fluid oz.)

Made in U.S.A.

Control Number on Bottom

Manufactured by:



**METREX RESEARCH  
CORPORATION**

Perle, Colorado 80134

ACCEPTED  
with COMMENTS  
in EPA Letter Dated:

NOV 10 1988

Under the Federal Insecticide,  
Fungicide, and Rodenticide Act,  
amended, for the pesticide  
registered under EPA Reg. No.

46781-5

**BEST AVAILABLE COPY**