



UNITED STATES ENVIRONMENTAL PROTECTION AGENCY
WASHINGTON, DC 20460

OFFICE OF CHEMICAL SAFETY
AND POLLUTION PREVENTION

June 16, 2023

Annette Marine
Agent
Silversan AG
Electronic Transmittal: annette@wagnerreg.com

Subject: Notification per PRN 98-10 – Update company name and address and general formatting.
Product Name: Aqua-Tec™
EPA Registration Number: 91025-1
Received Date: December 2, 2022
Action Case Number: 00414082

Dear Annette Marine:

The Agency is in receipt of your Application for Pesticide Notification under Pesticide Registration Notice (PRN) 98-10 for the above referenced product. The Antimicrobials Division (AD) has conducted a review of this request for its applicability under PRN 98-10 and finds that the action requested falls within the scope of PRN 98-10.

The label submitted with the application has been stamped “Notification” and will be placed in our records.

Should you wish to add/retain a reference to the company’s website on your label, then please be aware that the website becomes labeling under the Federal Insecticide Fungicide and Rodenticide Act and is subject to review by the Agency. If the website is false or misleading, the product would be misbranded and unlawful to sell or distribute under FIFRA section 12(a)(1)(E). 40 CFR 156.10(a)(5) lists examples of statements EPA may consider false or misleading. In addition, regardless of whether a website is referenced on your product’s label, claims made on the website may not substantially differ from those claims approved through the registration process. Therefore, should the Agency find or if it is brought to our attention that a website contains false or misleading statements or claims substantially differing from the EPA approved registration, the website will be referred to the EPA’s Office of Enforcement and Assurance.

If you have any questions, please contact Zebora Johnson at (202) 566-0730 or by mail at johnson.zebora@epa.gov

Sincerely,

A handwritten signature in cursive script that reads "Zebora Johnson".

For
Steven Snyderman
Product Manager, 33
Regulatory Management Branch II
Antimicrobials Division 97510M)

{MASTER LABEL}

{Front Panel}

Protects your lab samples
With confidence

Swizz Style, Inc.

Aqua-Tec™

Antimicrobial Water Preservation Cell for Use in CO2 Incubators and Water Baths
To Prevent Mold, Mildew, Algae, and Non-Public Health Bacteria

Laboratory Use Only

NOTIFICATION

91025-1

PREVENTS water contamination

The applicant has certified that no changes, other than those reported to the Agency have been made to the labeling. The Agency acknowledges this notification by letter dated:

Protects up to 6 months
Designed for all laboratory water types

06/16/2023

Active Ingredient:

Silver (as elemental) 62%

Other Ingredients: 38%

Total: 100%

Precautionary Statements and First Aid

See Instructions on Reverse

Keep Out of ~~the~~ Reach of Children
CAUTION

Net weight: _____

{Back Panel}

FIRST AID

| | |
|---------------------------------------|---|
| IF IN EYES | <ul style="list-style-type: none"> • Hold eye open and rinse slowly and gently with water for 15-20 minutes. • Remove contact lenses, if present, after the first 5 minutes, then continue rinsing eye. • Call poison control center or doctor for treatment advice. |
| IF SWALLOWED | <ul style="list-style-type: none"> • Immediately cCall a poison control center or doctor <u>immediately for treatment advice.</u> • <u>Have person sip a glass of water if able to swallow.</u> • Do not induce vomiting unless directed by a poison control center or doctor. • Do not give any liquid to the person. • Do not give anything by mouth to an unconscious person. |
| IF ON SKIN OR CLOTHING OR SKIN | <ul style="list-style-type: none"> • Take off contaminated clothing. • Rinse skin immediately with plenty of water for 15-20 minutes. • Call a poison control center or doctor for treatment advice. |
| IF INHALED | <ul style="list-style-type: none"> • <u>Move person to fresh air.</u> • <u>If person is not breathing, call 911 or an ambulance, then give artificial respiration, preferably by mouth-to-mouth, if possible.</u> |

| | |
|--|---|
| | <ul style="list-style-type: none"> • <u>Call a poison control center or doctor for treatment advice.</u> <p><u>Have the product container or label with you when calling a poison control center or doctor, or going to treatment.</u></p> <p><u>For medical emergencies call the poison control center at 1-800-222-1222.</u></p> |
|--|---|

FOR MEDICAL ASSISTANCE IN THE US, CALL 1-800-222-1222

PRECAUTIONARY STATEMENTS

Hazards to Humans and Animals

CAUTION. Causes eye irritation. May be harmful if swallowed. Avoid contact with skin, eyes or clothing. Wash thoroughly with soap and water after handling. Remove contaminated clothing and wash before reuse.

DIRECTIONS FOR USE:

It is a violation of Federal Law to use this product in a manner inconsistent with its labeling.

Directions for Use:

1. Place Aqua-Tec in water reservoir (pan or external bottle) of the CO2 incubator or water bath.
2. Use the suction cups to keep ~~the~~ Aqua-Tec in a fixed position.
3. Replace Aqua-Tec every 6 months to maintain effective protection against contamination. Use the peel off sticker as a reminder.

STORAGE AND DISPOSAL

STORAGE: Do not store together with: Strong oxidants, strong acids (nitric acid, sulphuric acid), strong lye, halogens, halogen connections, or concentrated hydrogen peroxide solutions.

DISPOSAL: Dispose of by wrapping the exhausted cartridge in newspaper and place in a trash can. Ensure compliance with all local and state regulations.

[Manufactured for:]

SILVERSAN AG Swizz Style, Inc.
Industriestrasse 21465 W. 3rd St.

CH-9400, Rorschach, Switzerland
Dover, OH 44622

EPA Reg. No. ~~88804-R~~91025-1

EPA Est. No. ~~88804-XX-XXX~~_____

1-877-663-7895

Catalog No. _____

Publication No. _____

Rev. _____